



Design of Experiments for optimal production of malaria vaccines with *Pichia pastoris*



Application
Note

#09

#01

#02

#03

#04

J. Fricke¹, K. Pohlmann¹, F. Tatge¹, R. Lang¹,
A. Ellert², B. Joksch², B. W. Faber³,
R. Luttmann¹

¹ HAW – Hamburg University of Applied
Sciences, Germany

² Sartorius Stedim Biotech GmbH,
Germany

³ BPRC – Biomedical Primate Research
Centre, The Netherlands

Application background

The aim of a joint project of the HAW Hamburg and the Biomedical Primate Research Center (BPRC) was to identify optimal protein expression conditions for high yields of potential malaria vaccines in the methylotrophic yeast *Pichia pastoris*.

Optimal expression conditions vary in accordance to the kind of *Pichia pastoris* strain used for the foreign protein expressed. As strain characterization and target protein expression optimization need a large number of cultivation runs, a fully automated multi-bioreactor plant was designed and built up for the implementation of investigations via Design of Experiments (DoE).


Enabling technology for future process development

Figure 1 shows the fully automated multi-bioreactor plant used for the executed DoE investigations. The plant design enables a reproducible inoculation of the six-fold parallel parameter screening reactor system BIOSTAT® Qplus with cell broth from the pre-cultivation bioreactor BIOSTAT® Bplus.

The process is operated with the SCADA system BioPAT® MFCS/win. Recipe structures following ANSI | ISA-88.01 permit automated cultivations including harvest, refresh and inoculation operations. The plant concept envisages DoE procedures, planned and carried out by the developed MFCS/win DoE module networked with the DoE software MODDE® by MKS Instruments AB.



Fig. 1: Multi-bioreactor DoE-plant with BIOSTAT® Bplus and Qplus

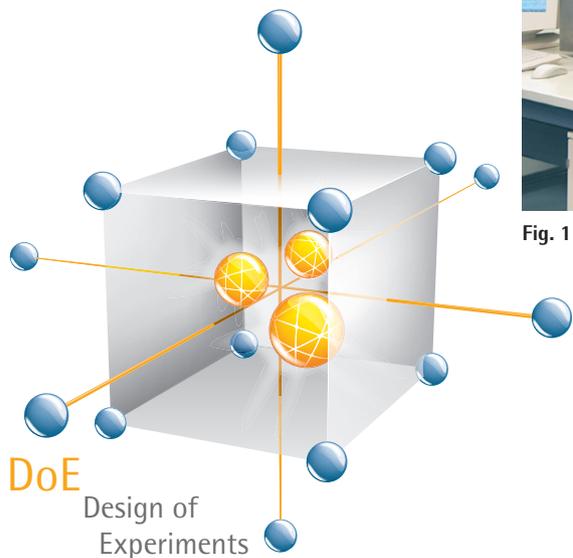

Investigated expression factors and quality responses

As a first step a DoE screening procedure focused on the impact of three parameters, the methanol concentration in the media phase c_{S2M} , the pH-value, and the liquid temperature ϑ_L during expression. By varying these process parameters with DoE it is expected to improve target protein yield with minimum proteolytic degradation.

Figure 2 shows the optimization potential of the chosen screening parameters by comparing two experiments. Experiment 1 with low pH-value and low expression temperature ϑ_L and a higher c_{S2M} level has a much lower target protein production c_{P1M} than experiment 2 with higher pH-value and ϑ_L and a lower c_{S2M} .


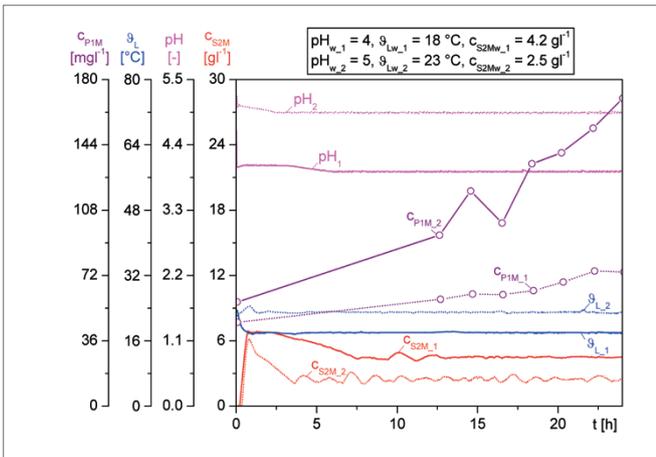


Fig. 2: Time course of two *Pichia pastoris* cultivations with different operation parameters reveals optimization potential

Creating a seamless experimental workflow

The wizard functionality of the MFCS/win DoE module was used to generate a set of screening experiments which were then automatically transferred into the already existing S88 recipe structure. The designed experiments were started simultaneously from within the DoE module and once triggered ran like a standard MFCS/win batch.

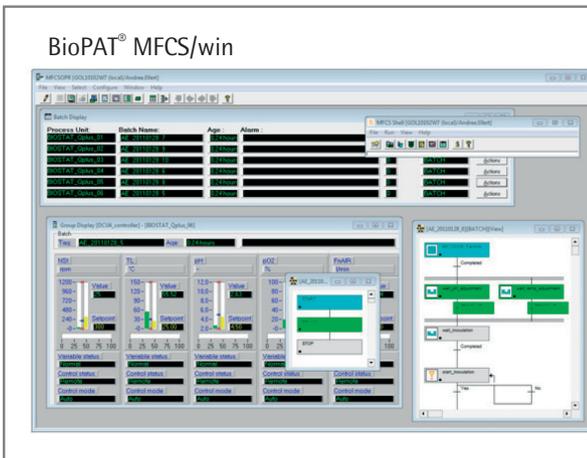

Results of the DoE-screening

Figure 4 shows the results of the DoE-screening which was accomplished by using a full factorial design. The lower response surface represents the productivity at pH 4, the upper at pH 6. The response surface plot indicates a higher productivity in the upper right corner with a low methanol concentration c_{S2M} , a higher temperature ϑ_L and a high pH-value.

Based on the screening results a more thorough DoE-procedure with optimization and robustness analyses are currently in progress. In the upper right corner of Fig. 4 the proposed CCC optimization is already shown.

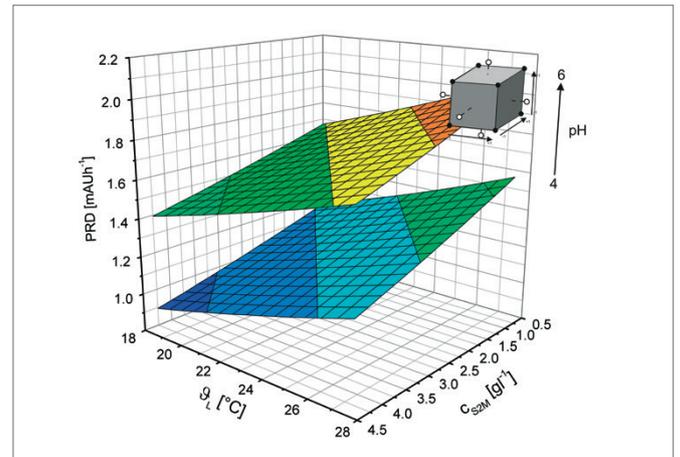


Fig. 4: Response surface plot of DoE-Screening of productivity (PRD) as a function of methanol concentration (c_{S2M}) and cultivation temperature (ϑ_L)

Following completion of the process, the DoE module extracts and stores the specified response data and automatically puts them into a MODDE® worksheet for further analysis, so that no typing errors and hence no odd anomalies hinder subsequent statistical data analysis.

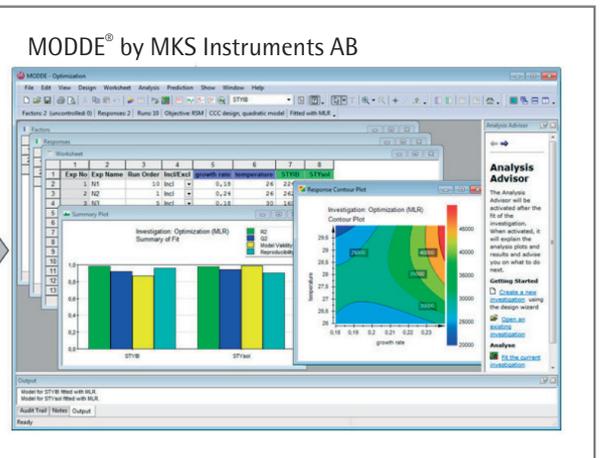


Fig. 3: Integration of DoE into the bioprocess management software BioPAT® MFCS/win

Sales and Service Contacts

For further contacts, visit www.sartorius-stedim.com

Europe

Germany

Sartorius Stedim Biotech GmbH
August-Spindler-Strasse 11
37079 Goettingen

Phone +49.551.308.0
Fax +49.551.308.3289

Sartorius Stedim Systems GmbH
Robert-Bosch-Strasse 5 – 7
34302 Guxhagen

Phone +49.5665.407.0
Fax +49.5665.407.2200

France

Sartorius Stedim FMT S.A.S.
ZI des Paluds
Avenue de Jouques – CS 91051
13781 Aubagne Cedex

Phone +33.442.845600
Fax +33.442.845619

Sartorius Stedim France SAS
ZI des Paluds
Avenue de Jouques – CS 71058
13781 Aubagne Cedex

Phone +33.442.845600
Fax +33.442.846545

Austria

Sartorius Stedim Austria GmbH
Modcenterstrasse 22
1030 Vienna

Phone +43.1.7965763.18
Fax +43.1.796576344

Belgium

Sartorius Stedim Belgium N.V.
Rue Colonel Bourg 105
1030 Bruxelles

Phone +32.2.756.06.80
Fax +32.2.756.06.81

Hungary

Sartorius Stedim Hungária Kft.
Kagyló u. 5
2092 Budakeszi

Phone +36.23.457.227
Fax +36.23.457.147

Italy

Sartorius Stedim Italy S.r.l.
Via dell'Antella, 76/A
50012 Antella-Bagno a Ripoli (FI)

Phone +39.055.63.40.41
Fax +39.055.63.40.526

Netherlands

Sartorius Stedim Netherlands B.V.

Phone +31.30.60.25.080
Fax +31.30.60.25.099

filtratie.nederland@sartorius-stedim.com

Poland

Sartorius Stedim Poland Sp. z o.o.
ul. Wrzesinska 70
62-025 Kostrzyn

Phone +48.61.647.38.40
Fax +48.61.879.25.04

Russian Federation

LLC "Sartorius Stedim RUS"
Vasilyevsky Island
5th line 70, Lit. A
199178 St. Petersburg

Phone +7.812.327.53.27
Fax +7.812.327.53.23

Spain

Sartorius Stedim Spain, S.A.U.
Avda. de la Industria, 32
Edificio PAYMA
28108 Alcobendas (Madrid)

Phone +34.913.586.098
Fax +34.913.589.623

Switzerland

Sartorius Stedim Switzerland AG
Ringstrasse 24 a
8317 Tagelswangen

Phone +41.52.354.36.36
Fax +41.52.354.36.46

U.K.

Sartorius Stedim UK Ltd.
Longmead Business Centre
Blenheim Road, Epsom
Surrey KT19 9 QQ

Phone +44.1372.737159
Fax +44.1372.726171

Ukraine

LLC "Sartorius Stedim RUS"
Post Box 440 "B"
01001 Kiev, Ukraine

Phone +380.44.411.4918
Fax +380.50.623.3162

Americas

USA

Sartorius Stedim North America Inc.
5 Orville Drive, Suite 200
Bohemia, NY 11716

Toll-Free +1.800.368.7178
Fax +1.631.254.4253

Argentina

Sartorius Argentina S.A.
Int. A. Ávalos 4251
B1605ECS Munro
Buenos Aires

Phone +54.11.4721.0505
Fax +54.11.4762.2333

Brazil

Sartorius do Brasil Ltda
Avenida Senador Vergueiro 2962
São Bernardo do Campo
CEP 09600-000 – SP- Brasil

Phone +55.11.4362.8900
Fax +55.11.4362.8901

Mexico

Sartorius de México, S.A. de C.V.
Libramiento Norte de Tepotzotlan s/n,
Colonia Barrio Tlacateco,
Municipio de Tepotzotlan,
Estado de México,
C.P. 54605

Phone +52.55.5562.1102
Fax +52.55.5562.2942

leadsmex@sartorius.com

Peru

Sartorius Peru S.A.C.
Avenue Alberto del Campo 411
Floor 12 – The Office
15076 – San Isidro, Lima

Phone +51.1.441 0158
Fax +51.1.422 6100

Asia | Pacific

Australia

Sartorius Stedim Australia Pty. Ltd.
Unit 5, 7-11 Rodeo Drive
Dandenong South Vic 3175

Phone +61.3.8762.1800
Fax +61.3.8762.1828

China

Sartorius Stedim Biotech (Beijing) Co. Ltd.
No. 33 Yu'an Road
Airport Industrial Park Zone B
Shunyi District, Beijing 101300

Phone +86.10.80426516
Fax +86.10.80426580

Sartorius Stedim (Shanghai)
Trading Co., Ltd.
3rd Floor, North Wing, Tower 1
No. 4560 Jinke Road
Zhangjiang Hi-Tech Park
Pudong District
Shanghai 201210, P.R. China

Phone +86.21.6878.2300
Fax +86.21.6878.2882

Sartorius Stedim Biotech (Beijing) Co. Ltd.
Guangzhou Representative Office
Unit K, Building 23
Huihua Commerce & Trade Building
No. 80 Xianlie Middle Road
Guangzhou 510070

Phone +86.20.37618687 | 37618651
Fax +86.20.37619051

India

Sartorius Stedim India Pvt. Ltd.
#69/2-69/3, NH 48, Jakkasandra
Nelamangala Tq
562 123 Bangalore, India

Phone +91.80.4350.5250
Fax +91.80.4350.5253

Japan

Sartorius Stedim Japan K.K.
4th Fl., Daiwa Shinagawa North Bldg.
8-11, Kita-Shinagawa 1-chome
Shinagawa-ku, Tokyo, 140-0001 Japan

Phone +81.3.4331.4300
Fax +81.3.4331.4301

Malaysia

Sartorius Stedim Malaysia Sdn. Bhd.
Lot L3-E-3B, Enterprise 4
Technology Park Malaysia
Bukit Jalil
57000 Kuala Lumpur, Malaysia

Phone +60.3.8996.0622
Fax +60.3.8996.0755

Singapore

Sartorius Stedim Singapore Pte. Ltd.
10 Science Park Rd
The Alpha #02-13/14
Singapore Science Park II
Singapore 117684

Phone +65.6872.3966
Fax +65.6778.2494

South Korea

Sartorius Korea Biotech Co., Ltd.
8th Floor, Solid Space B/D,
PanGyoYeok-Ro 220, BunDang-Gu
SeongNam-Si, GyeongGi-Do, 463-400

Phone +82.31.622.5700
Fax +82.31.622.5799



▶ www.sartorius-stedim.com